


National Diagnostic Protocol

Potato Mop-Top Virus (PMTV)

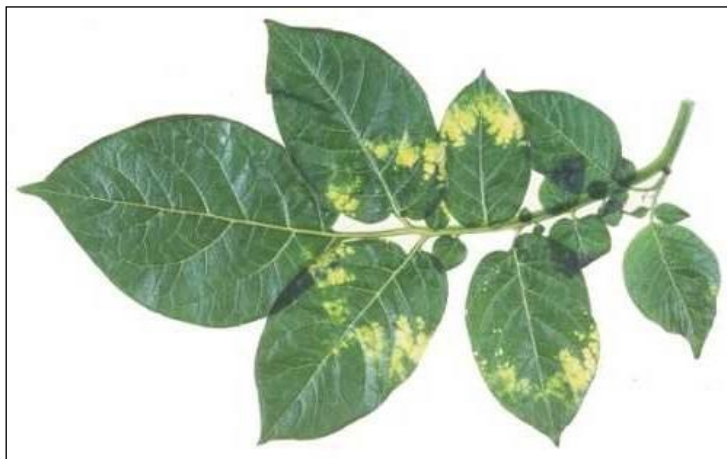

NDP 15 V2

© Commonwealth of Australia

Ownership of intellectual property rights

Unless otherwise noted, copyright (and any other intellectual property rights, if any) in this publication is owned by the Commonwealth of Australia (referred to as the Commonwealth).

Creative Commons licence

All material in this publication is licensed under a Creative Commons Attribution 3.0 Australia Licence, save for content supplied by third parties, logos and the Commonwealth Coat of Arms.



Creative Commons Attribution 3.0 Australia Licence is a standard form licence agreement that allows you to copy, distribute, transmit and adapt this publication provided you attribute the work. A summary of the licence terms is available from <http://creativecommons.org/licenses/by/3.0/au/deed.en>. The full licence terms are available from <https://creativecommons.org/licenses/by/3.0/au/legalcode>.

This publication (and any material sourced from it) should be attributed as: Subcommittee on Plant Health Diagnostics (2020). National Diagnostic Protocol for Potato Mop-Top Virus (PMTV) – NDP15 V2. (Eds. Subcommittee on Plant Health Diagnostics) Authors Eyre J., Rowles-van Rijswijk B., Rodoni B. and Lovelock D.; Reviewers Thomas J. and Lovelock D. ISBN 978-0-6481143-7-6 CC BY 3.0.

Cataloguing data

Subcommittee on Plant Health Diagnostics (2020). National Diagnostic Protocol for Potato Mop-Top Virus (PMTV) - NDP15 V2. (Eds. Subcommittee on Plant Health Diagnostics) Authors Eyre J., Rowles-van Rijswijk B., Rodoni B. and Lovelock D.; Reviewers Thomas J. and Lovelock D. ISBN 978-0-6481143-7-6.

ISBN; 978-0-6481143-7-6.

Internet

Report title is available at: <http://plantbiosecuritydiagnostics.net.au/resource-hub/priority-pest-diagnostic-resources/>

Department of Agriculture and Water Resources

Street Address 1: [18 Marcus Clarke Street](#), Canberra City ACT 2601

Street Address 2: [7 London Circuit](#), Canberra City ACT 2601

Postal Address: GPO Box 858, Canberra City ACT 2601

Switchboard Phone: 02 6272 3933

Web: <http://www.agriculture.gov.au>

Inquiries regarding the licence and any use of this document should be sent to: copyright@agriculture.gov.au.

The Australian Government acting through the Department of Agriculture and Water Resources has exercised due care and skill in the preparation and compilation of the information and data in this publication. Notwithstanding, the Department of Agriculture and Water Resources, its employees and advisers disclaim all liability, including liability for negligence, for any loss, damage, injury, expense or cost incurred by any person as a result of accessing, using or relying upon any of the information or data in this publication to the maximum extent permitted by law.

Purpose

National Diagnostic Protocols (NDPs) are diagnostic protocols for the unambiguous taxonomic identification of plant pests. NDPs:

- are a verified information resource for plant health diagnosticians
- are consistent with ISPM No. 27 – Diagnostic Protocols for Regulated Pests
- provide a nationally consistent approach to the identification of plant pests enabling transparency when comparing diagnostic results between laboratories; and,
- are endorsed by regulatory jurisdictions for use (either within their own facilities or when commissioning from others) in a pest incursion.

Where an International Plant Protection Convention (IPPC) diagnostic protocol exists it should be used in preference to NDPs although NDPs may contain additional information to aid diagnosis. IPPC protocols are available on the IPPC website:

<https://www.ippc.int/core-activities/standards-setting/ispms>

Process

NDPs are facilitated and endorsed by the Subcommittee on Plant Health Diagnostics (SPHD). SPHD reports to Plant Health Committee and is Australia's peak technical and policy forum for plant health diagnostics.

NDPs are developed and endorsed according to Reference Standards developed and maintained by SPHD. Current Reference Standards are available at

<http://plantbiosecuritydiagnostics.net.au/sphd/sphd-reference-standards/>

NDPs are living documents. They are updated every 5 years or before this time if required (i.e. when new techniques become available).

Document status

This version of the National Diagnostic Protocol (NDP) for Potato Mop-Top Virus (PMTV) is current as at the date contained in the version control box below.

| | |
|-----------------|--------------------------|
| PEST STATUS | Not present in Australia |
| PROTOCOL NUMBER | NDP 15 |
| VERSION NUMBER | V2 |
| PROTOCOL STATUS | Endorsed |
| ISSUE DATE | 2020 |
| REVIEW DATE | 2025 |
| ISSUED BY | SPHD |

The most current version of this document is available from the SPHD website:

<http://plantbiosecuritydiagnostics.net.au/resource-hub/priority-pest-diagnostic-resources/>

Further information

Inquiries regarding technical matters relating to this project should be sent to:

sphds@agriculture.gov.au

Contents

| | | |
|----------|--|-----------|
| 1 | INTRODUCTION..... | 2 |
| 1.1 | Host range..... | 2 |
| 1.2 | Transmission | 2 |
| 2 | TAXONOMIC INFORMATION..... | 3 |
| 3 | DETECTION..... | 4 |
| 3.1 | Diagnostic flow chart..... | 4 |
| 3.2 | Symptom description..... | 5 |
| 3.3 | Diseases causing symptoms similar to <i>Potato mop-top virus</i> | 9 |
| 3.4 | Sample Collection..... | 10 |
| 4 | IDENTIFICATION..... | 11 |
| 4.1 | Molecular methods..... | 11 |
| 4.2 | Serological Tests – ELISA..... | 19 |
| 5 | CONTACTS FOR FURTHER INFORMATION..... | 20 |
| 6 | ACKNOWLEDGEMENTS..... | 21 |
| 7 | REFERENCES..... | 22 |
| 7.1 | Related Articles..... | 24 |
| 8 | APPENDIX..... | 26 |
| 8.1 | ELISA Buffers | 26 |
| 9 | DIAGNOSTICS PROCEDURES TO SUPPORT SURVEILLANCE | 27 |
| 9.1 | Introduction | 27 |
| 9.2 | Sampling..... | 27 |
| 9.3 | In Field Tests..... | 27 |
| 9.4 | Laboratory Tests | 27 |
| 9.5 | Acknowledgements..... | 28 |
| 9.6 | References..... | 28 |

1 INTRODUCTION

PMTV produces internal and external potato tuber symptoms, seriously affecting tuber quality and reducing yield (Germundsson *et al.*, 2002). Infection with PMTV causes some quantitative yield loss but the qualitative losses are more important and can lead to total crop rejection by supermarkets and processors (Jones and Harrison, 1972; Mumford *et al.*, 2000; Yellareddygaru *et al.*, 2018).

1.1 Host range

PMTV affects potato (*Solanum tuberosum*).

Weeds from the Chenopodiaceae and Solanaceae are susceptible

(<https://link.springer.com/article/10.1007/BF02732217>, Anderson *et al.* 2002), Anderson *et al.* 2002).

1.2 Transmission

The plasmodiophoromycete *Spongospora subterranea* f. sp. *subterranea* is the only known vector of PMTV (Helias *et al.*, 2003; Sokmen *et al.*, 1998). *S. subterranea* transmits the virus to potato roots via viruliferous fungal zoospores (Sokmen *et al.*, 1998). The vector of PMTV is also the causal agent of powdery scab, and has been associated with potato production since the mid-1800s (Büchen-Osmond, 2001 onwards). PMTV can persist in the soil for many years in the resting spores of its vector (Helias *et al.*, 2003; Santal *et al.*, 2010), and soils therefore can remain infective for extended periods of time (Sokmen *et al.*, 1998). It has been shown that PMTV can persist for at least two years in the resting spores of *S. subterranea* and, PMTV was also detected in field soil 12 years after the last potato crop had been grown in that soil (Jones and Harrison, 1972). Soil type does not influence the occurrence of PMTV, as the virus has been detected in soil types ranging from light sands to heavy loams (Jones and Harrison, 1972).

A number of plant families are susceptible to infection by powdery scab. Under greenhouse conditions plant species within the *Solanaceae*, *Chenopodiaceae* and *Cruciferae* were found to be hosts of *S. subterranea* (Jones and Harrison, 1972). However, Johnson (2004) reported that only potato and black nightshade (*Solanum nigrum*) produced long-lived fungal resting spores of *S. subterranea*. More recently, Tsror *et al.*, (2020) observed *S. subterranean* successfully infecting a wider weed host range including species within the families of *Malvaceae* and *Zygophyllaceae*.

2 TAXONOMIC INFORMATION

Species *Potato mop top virus*

Genus *Pomovirus*

Family *Virgaviridae*

Order *Martellivirales*

Class *Alsuviricetes*

Phylum *Kitrinoviricota*

Kingdom *Orthornavirae*

Realm *Riboviria*

PMTV is the type member of the *Pomovirus* genus (Büchen-Osmond, 2001 onwards; Cerovska *et al.*, 2003; Tidona and Darai, 2002), and has straight, tubular, rigid particles, measuring 18-20 nm in diameter and 100-300 nm in length (Figure) (Cerovska *et al.*, 2003; Tidona and Darai, 2002). Discrepancies in the reported length of the viral particle are most probably due to the fragility of the particles, which readily disintegrate (Cerovska *et al.*, 2003). PMTV occurs in the hosts' cytoplasm (Tidona and Darai, 2002), and is unevenly distributed and present at low concentration in infected tissues (Helias *et al.*, 2003).

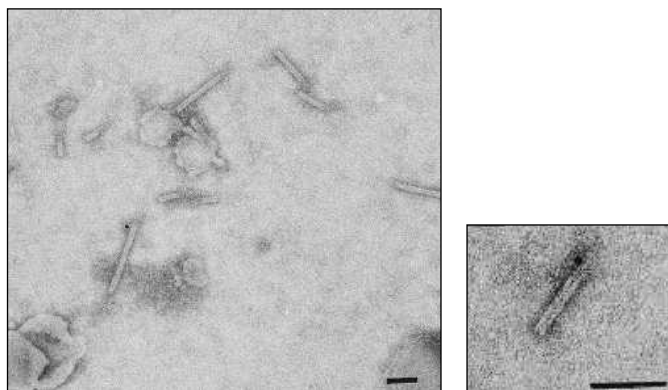



Figure 1. PMTV particles (left), uncoiling of the PMTV particle (right). In each electron micrograph, the bar represents 100 nm (Cowan *et al.*, 1997).

The PMTV genome consists of three positive-sense, single-stranded RNA molecules; RNA1 (6 kb), RNA2 (3.2 kb) and RNA3 (2.5 kb) (Helias *et al.*, 2003; Mayo *et al.*, 1996; Sokmen *et al.*, 1998; Tidona and Darai, 2002). RNA 1 encodes a methyltransferase, a helicase, and the viral RNA dependent RNA polymerase. RNA 2 encodes the triple gene block and an 8K, cysteine-rich protein of unknown function. RNA 3 encodes coat protein and readthrough protein (Nielsen and Nicolaisen, 2003). Comparisons of the PMTV coat protein nucleotide sequence of various isolates has shown that the coat protein open reading frame is highly conserved (Helias *et al.*, 2003). The 3'- and 5'-terminal sequences of PMTV are highly conserved and could prove useful in the design of specific primers (Kalyandurg *et al.*, 2017; Gil *et al.*, 2016).

3 DETECTION

Please note, that the PMTV detection system outlined in this manual is based on the RT-PCR protocols published by Sokmen *et al.* (1998) and Mayo *et al.* (1996). An additional nested PCR (Zhou *et al.*, 2019) can also be used.

3.1 Diagnostic flow chart

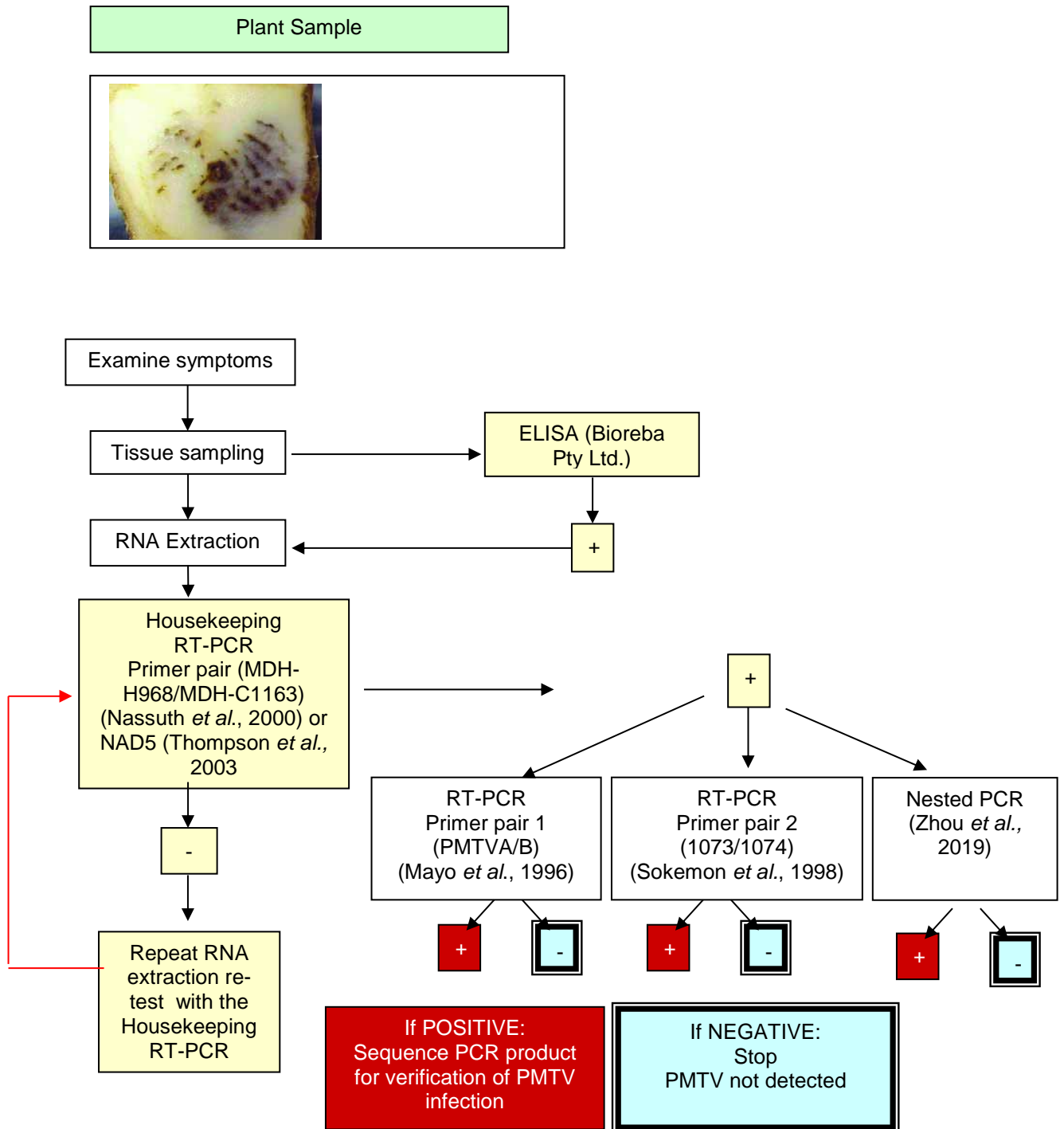

3.2 Symptom description

PMTV symptoms on potato include ‘mop-top’ stunting of foliage (Figure), yellow blotching (Figure) and yellow V-shaped markings (**Error! Reference source not found.**) on potato leaves, and sporadic shortening of internodes (Figure 5) (Calvert, 1968; Mayo *et al.*, 1996; Scott *et al.*, 1994).

PMTV causes a wide range of symptoms in potato haulms (stems) and tubers depending on viral strain, environmental conditions and cultivar susceptibility (Kurppa, 1989). Studies have shown that PMTV symptom expression can vary in the same potato variety, and several virus isolates differing in virulence have been reported (Harrison and Jones, 1970; Mayo *et al.*, 1996; Nielsen and Nicolaisen, 2003; Santala *et al.*, 2010; Gil *et al.*, 2016).

Tubers infected with PMTV often display rust coloured, necrotic arcs, flecks, rings or lines through the tuber flesh (Figure , 11, 12, 13). The brown necroses often penetrate into the tuber’s flesh almost without interruption (**Error! Reference source not found.**). These symptoms are known as “spraing”, or “corky ring-spot”. Spraing symptoms (brown lines and arcs in the tuber flesh) in susceptible cultivars are more prevalent in cool weather (around 15°C) (Sokmen *et al.*, 1998). External symptoms include cracking (Figure), raised necrotic rings (Figure) and raised necrotic lines (**Error! Reference source not found.**). Tuber symptoms may be absent at harvest, but can develop during storage. Visual diagnosis based on spraing induced by PMTV is difficult as *Tobacco rattle virus* (TRV), *Potato virus Y NTN* strain (PVY^{NTN}) and internal rust spot can also induce similar symptoms (Figures 14-17).



Figure 2. Healthy potato haulm on the left and PMTV infected potato on right showing “mop top” symptoms (© Scottish Agricultural Science Agency).



Figure 3. Yellow blotching on leaves (© Scottish Agricultural Science Agency).

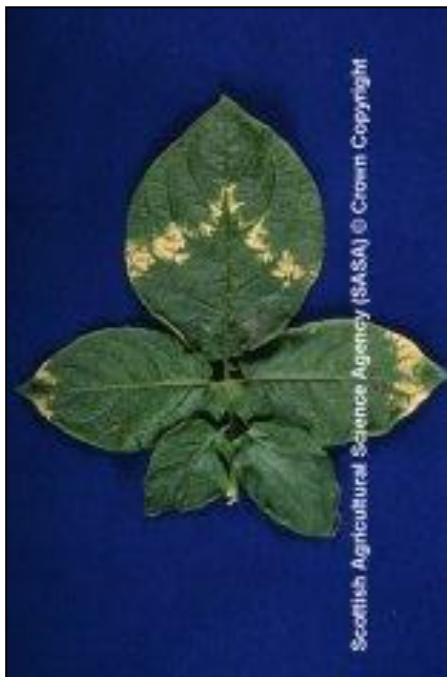


Figure 4. Chlorotic V-shaped markings in the leaflets (© Scottish Agricultural Science Agency)



Figure 5. Sporadic shortening of internodes (source: <http://www.unece.org/trade/agr/standard/potatoes/PestPicturesE.htm>).

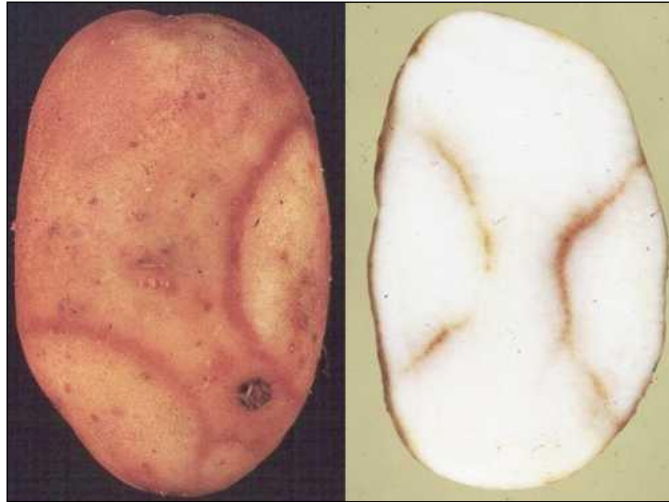


Figure 6. External (left) and internal (right) PMTV spraing symptoms in naturally infected potato tubers cv. Arran Pilot (© Scottish Agricultural Science Agency).



Figure 7. Cracked and distorted naturally infected tubers of potato cv. Alpha produced in second year of infection (© Scottish Agricultural Science Agency).



Figure 8. Tuber cracking and scaring (© Scottish Agricultural Science Agency).



Figure 9. Raised necrotic rings on tuber (© Scottish Agricultural Science Agency).



Figure 10. Raised necrotic lines (© Scottish Agricultural Science Agency).



Figure 11. Multiple necrotic concentric arcs in tuber flesh (© Johnson, University of Maine).



Figure 12. Necrotic arcs in tuber flesh (© Johnson, University of Maine).



Figure 13. Necrotic arcs on potato crisps, UK (Source: Rodoni, DPI Victoria).

3.3 Diseases causing symptoms similar to *Potato mop-top virus*

There are several viruses which produce symptoms similar to PMTV. *Tobacco rattle virus* (TRV) can cause yellow blotches on leaves () and necrotic arcs in tubers () similar to PMTV. The tuber necrosis strain of *Potato virus Y* (PVY^{NTN}) causes necrotic rings to appear on potato tubers (**Error! Reference source not found.**). The physiological disorder, “internal rust spot” can also cause tuber symptoms (**Error! Reference source not found.**) easily confused with PMTV.



Figure 14. Potato showing leaf yellowing symptoms of infection by TRV (Source: British potato council).

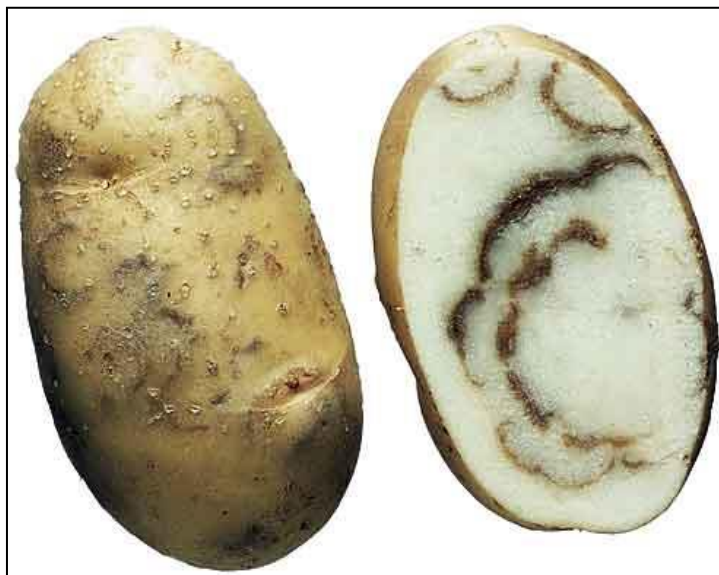


Figure 1. Potato tuber showing necrotic arcs in the flesh caused by infection by TRV (Source: British potato council, http://www.potato.org.uk/department/knowledge_transfer/pests_and_diseases/ref.html?item=27.)

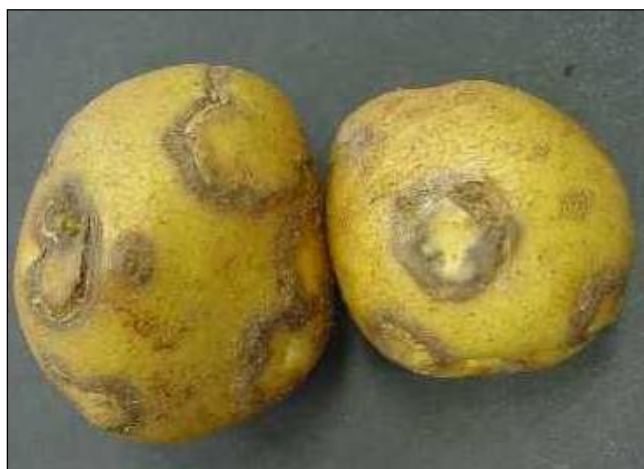


Figure 2. Tubers infected with Potato Tuber Necrotic Ring Spot Disease (PVY^{NTN}) (Source: Dr Brendan Rodoni, DPI-Victoria).



Figure 17. Tuber symptoms caused by internal rust spot (Source: British potato council).

3.4 Sample Collection

Leaf and tuber samples should not have free water on their surfaces. Samples should not be collected during hot weather and all samples should be cooled and transported to a testing facility in a cool environment if possible.

3.4.1 Leaf sampling

Record plant symptoms. Older symptomatic leaves around the middle of the haulm (stem) (Figure , 6), and leaves showing unusual symptoms must be sampled. Store plant material at 4°C or at -20°C until required.

3.4.2 Tuber sampling

Tubers removed from cold storage should be incubated at 20°C for four weeks prior to testing. This incubation has shown to double the sensitivity of detection (Sokmen *et al.*, 1998). If fresh tubers are used negative test results should be treated with caution due to the low levels of virus in the tuber. Stolon and rose ends of the tuber (Figure) must be included in each sample. PMTV does not infect all the tubers produced by an infected mother plant and therefore only a variable proportion of the progeny plants generated from the tuber seed are infected with PMTV in the following year (Sokmen *et al.*, 1998).

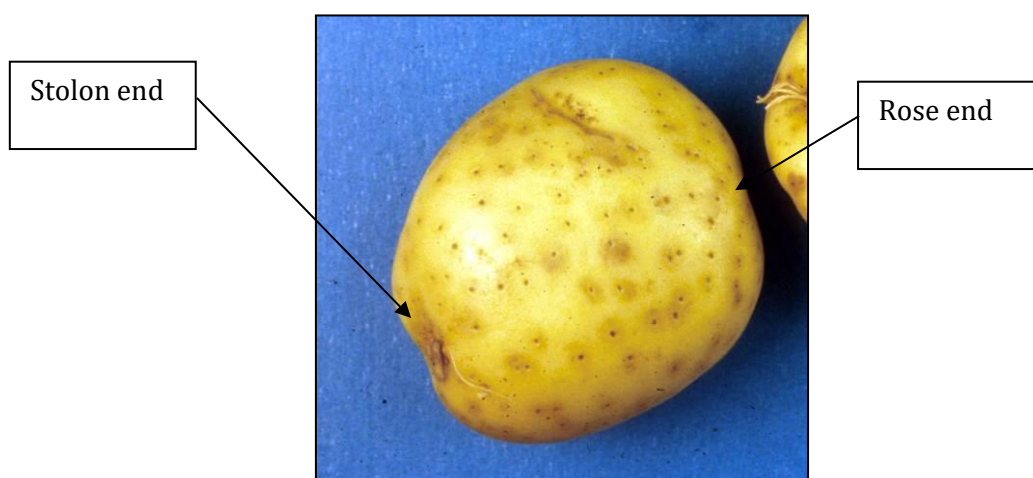


Figure 183. Stolon and rose end of tuber (Source: B Hall, SARDI)

...

4 IDENTIFICATION

4.1 Molecular methods

4.1.1 *Total RNA extraction*

Wear disposable gloves and a lab coat at all times. Have an autoclave bag ready to dispose of all plant material, tips, tubes, gloves and paper towel that have come into contact with any suspect plant material. A footbath containing disinfectant located at the doorway of the laboratory must be used when exiting the lab.

Equipment required

1. 2-20 μL , 20-200 μL , and 200-1000 μL micropipettes and sterile tips
2. Autoclave
3. Autoclave bags
4. Balance (at least 2 decimal places)
5. Disposable gloves
6. Disposable plastic pasteur pipettes
7. Microcentrifuge
8. RNeasy® Plant Mini Kit (Qiagen™) (Or similar RNA extraction Kit)
9. Sterile microcentrifuge tubes
10. Paper towel
11. Sharps container
12. Sterile scalpel blades and scalpel blade handle
13. Waterbath or heatblock set at 70°C
14. Weighboats

ALSO

- 1a: Autoclaved mortar and pestle, Fume hood, sterile sand

OR

- 1b: Qiagen Tissue Lyser(Qiagen™), 2 ml snap-lock tubes, Stainless steel beads, Liquid Nitrogen

OR

- 1c: Homex tissue macerater, Homex bags (Bioreba AG / BioSys), Plastic disposable pasteur pipettes

Reagents required

1. MacKenzie buffer (MacKenzie *et al.*, 1997)

| Chemical | Amount | Final Concentration |
|--|---------|---------------------|
| Guanidine thiocyanate (CH ₅ N ₃ ·CHNS) | 23.64 g | 4 M |
| 3M Sodium acetate (C ₂ H ₃ NaO ₂) | 3.33 ml | 0.2 M |
| 0.5M EDTA (C ₁₀ H ₁₆ N ₂ O ₈) | 2.5 ml | 25 mM |
| PVP-40 (Polyvinylpyrrolidone) | 1.25 g | 2.5% (w/v) |

Add sterile distilled water to final volume of 50 ml

Store at room temperature

Please note, fresh MacKenzie buffer should be prepared every 3-6 months

2. β-mercaptoethanol (C₂H₆OS)
3. 20% **N-Lauroylsarcosine solution (w/v)**
4. 100% Ethanol

Method

The RNA extraction method is based on that described by MacKenzie *et al.* (1997). All steps are carried out at room temperature as follows:

1a. If using a mortar and pestle to homogenise samples:

- 1a-1. Determine the number of samples and label plastic tubes accordingly.
- 1a-2. Use new clean gloves and scalpel blades for each sample.
- 1a-3. Cut each new sample on fresh paper towel on the bench.
- 1a-4. Weigh out 400 mg of plant sample (leaf or tuber).
- 1a-5. Place sample in mortar.
- 1a-6. Add 1980 µl of MacKenzie buffer.
- 1a-7. Add 20 µl of β-mercaptoethanol in the fumehood.
- 1a-8. Homogenise in fume hood.
- 1a-9. Pipette 1.0 ml of the mixture into a labelled microcentrifuge tube (you may need to cut the end of the pipette tip if the slurry is too thick).
- 1a-10. Continue to step 2.

1b. If using the Qiagen™ Tissue Lyser:

- 1b-1. Determine the number of samples and label the 2 ml snap-lock tubes accordingly.
- 1b-2. Use new clean gloves and scalpel blades for each sample.
- 1b-3. Cut each new sample on fresh paper towel on the bench.
- 1b-4. Weigh out 100 mg of plant sample (leaf or tuber) and place sample in the appropriate tube.
- 1b-5. Add 990 µl of MacKenzie buffer.
- 1b-6. Add 10 µl of β-mercaptoethanol in the fumehood.
- 1b-7. Close tubes.
- 1b-8. Place tubes in the Adaptor Set, in the Qiagen™ Tissue Lyser. Homogenise for 1 min at 30 Hz.
- 1b-9. Rotate tubes within the Adaptor Set, so that tubes in the centre are moved to the outside.
- 1b-10. Continue to step 2.

1c. If using the Homex tissue macerater:

- 1c-1. Determine the number of samples and label plastic tubes accordingly.
- 1c-2. Use new clean gloves and scalpel blades for each sample.
- 1c-3. Cut each new sample on fresh paper towel on the bench.
- 1c-4. Weigh out 200 mg of plant sample (leaf or tuber).
- 1c-5. Place sample in Homex bag.
- 1c-6. Add 1980 µl of MacKenzie buffer.
- 1c-7. Add 20 µl of β-mercaptoethanol in the fumehood.
- 1c-8. Macerate tissue with the Homex.
- 1c-9. With a plastic disposable pasteur pipette, transfer 1.0 ml of the mixture into a labelled microcentrifuge tube.
- 1c-10. Continue to step 2.

2. Carefully read the RNeasy® Mini Handbook.
3. Add 100 µL of 20% Sarkosyl to each tube and mix.
4. Incubate tubes at 70°C for 10 minutes.
5. Spin tubes in microcentrifuge for 1 minute at 13,000 rpm.
6. Continue with step 4 of the “RNeasy® Mini Protocol for Isolation of Total RNA from Plant Cells and Tissues and Filamentous Fungi” in the RNeasy® Mini Handbook and follow as per manufacturer's instructions.

4.1.2 Detection of PMTV in total RNA extracts using one step RT-PCR

For the reliable detection of PMTV, total RNA extracts are subjected to four RT-PCR tests, as outlined below, with primer sequences and annealing temperatures listed in Table 1 and 2.

1. **Primer pair PMTVA (R) / PMTVB (R).** This primer pair will detect the pathogen by amplifying approximately 550bp of the PMTV coat protein gene on RNA 3 (Mayo *et al.*, 1996).
2. **Primer pair 1073 (F) / 1074 (R).** This primer pair anneals to RNA 2 of the PMTV genome and amplifies 283bp of the 3' end of the triple-gene-block and the 8K protein genes (Sokmen *et al.*, 1998).
3. **Nested PCR:** 1st Round Primer pair H360 (F) / C819 (R) and 2nd Round Primer pair PMTV-171 (F) / PMTV-433 (R). These primer pairs anneal to the PMTV coat protein, resulting in amplicons at sizes of 460bp and 263bp respectively (Zhou *et al.*, 2019)
4. **House-keeping gene (Primer pair MDH-H968 (F) / MDH-C1163 (R).** The MDH-H968/MDH-C1163 primer pair are designed to amplify a 196 bp region of the plant mRNA encoding malate dehydrogenase (MDH) gene (Nassuth *et al.*, 2000). This gene is highly conserved among plants and therefore RT-PCR amplification of the MDH mRNA is used as an internal RT-PCR control to, a) determine the quality of the RNA extract, and b) determine whether the RNA extract contains inhibitors that will interfere with the activity of the reverse transcriptase and *Taq* DNA polymerase enzymes. Alternatively, the Primer pair of **NAD2.1a / Nad 2.2b** can be used to detect mRNA of the NADH dehydrogenase ND2 subunit; the amplicon size is approximately 188bp (Thompson *et al.*, 2003).

Equipment required

1. 0-2 μ L, 2-20 μ L, 20-200 μ L, & 200-1000 μ L pipettes and sterile tips
2. 0.2 or 0.5 mL sterile PCR tubes
3. Bulb spinner or centrifuge
4. Disposable gloves
5. Freezer
6. Gel electrophoresis tanks, rigs and racks
7. DNA Molecular Weight markers
8. Crushed ice
9. Sterile microcentrifuge tubes to store reagents
10. Microwave
11. Power pack
12. Thermocycler
13. UV transilluminator with camera

Reagents

1. Primers

For the detection of PMTV three specific primers sets are required (Table 1). Each primer is used as a stock solution at a concentration of 10 μ M.

Table 1 Primers required for the detection of PMTV

| Primer | Sequence (5'-3') | Target | Anneal temp. | PCR product size |
|-------------------------------|--------------------------------------|---|--------------|------------------|
| PMTVA (F) ¹ | CTA TGC ACC AGC CCA GCG T | PMTV RNA-3, nucleotide 268 to 801 | 50°C | 533 bp |
| PMTVB (R) ¹ | TCT CGG ATA CCA CCC TT | | | |
| 1073 (F) ² | AGT AGC AAG TAC GCC CTG TG | PMTV RNA-2, nucleotide 2419 to 2702 | 50°C | 283 bp |
| 1074 (R) ² | CCG AAT TCC TGT AAG CAC TAA CAC | | | |
| MDH- H968(F) ³ | GCA TCT GTG GTT CTT GCA GG | mRNA encoding malate dehydrogenase (MDH) | 54°C | 196 bp |
| MDH-C1163 (R) ³ | CCT TTG AGT CCA CAA GCC AA | | | |
| Nad2.1a ⁴ | GGA CTC CTG ACG TAT ACG AAG GAT C | NADH dehydrogenase ND2 subunit | 50°C | 188bp |
| Nad2.2b ⁴ | AGC AAT GAG ATT CCC CAA TAT CAT | | | |

¹(Mayo *et al.*, 1996), ²(Sokmen *et al.*, 1998), ³(Nassuth *et al.*, 2000) ⁴(Thompson *et al.*, 2003).

Table 2 Primers required for the detection of PMTV using a nested-PCR

| Primer | Sequence (5'-3') | Target | Anneal temp. | PCR product size |
|--------------|-----------------------------------|--|--------------|------------------|
| H360 (F) | CAT GAA GGC TGC CGT GAG GAA GT | PMTV Coat Protein (1 st Round) | 58°C | 460bp |
| C819 (R) | CTA TGC ACC AGC CCA GCG TAA CC | | | |
| PMTV-171 (F) | TAG CGT GTT GAG CCA GGT GT | PMTV Coat Protein (2 nd Round) | 58°C | 263bp |
| PMTV-433 (R) | CCT CTG CGA GTT GAT GTG CC | | | |

(Zhou *et al.*, 2019)

2. PCR Controls

1. Positive control - RNA extract from plant tissue infected with PMTV.
- Alternatively a “plasmid control” that has the target PMTV sequence cloned into the plasmid. (Positive controls available from DSMZ or contact author for further information).
2. Negative plant control - RNA extract from uninfected plant tissue of the same species as that used for the positive control.
3. Negative buffer control - an aliquot of the RT-PCR “Master Mix” without template.
4. The house keeping RT-PCR, using primers MDH-H968/MDH-C1163 or Nad2.1a/Nad2.2b, reduces the risk of false negative results. The generation of a band confirms the presence of RNA in the extract, and that the RNA extract does not contain inhibitors. Failure to produce an amplicon of expected size (196/188bp bp respectively) indicates that either dilution of the RNA extract or re-extraction of RNA from the sample is required.

3. RT-PCR reagents

1. One-step RT-PCR kit/PCR Kit (Kits validated within laboratories can also be used):
 - Invitrogen® SuperScript™ III One-Step RT-PCR with Platinum® *Taq* DNA Polymerase, Catalogue No. 12574026, is recommended.
 - Qiagen One step RT –PCR kit (Cat. No. 210212) was successfully used in verification of protocol while following the manufacturer’s instructions.
2. For the nested PCR Invitrogen® *Taq* DNA Polymerase Catalogue No. 109660108 is recommended for the second PCR. Nuclease-free water

4. 5x TBE Buffer

| | Per 1 litre |
|---|-------------|
| Tris (C ₄ H ₁₁ NO ₃) | 54 g |
| Boric acid (H ₃ BO ₃) | 27.5 g |
| 0.5M EDTA ([CH ₂ N(CH ₂ .COOH).CH ₂ COONa] ₂ .2H ₂ O) pH 8.0 | 20 ml |

Store at room temperature.

5. SYBR safe (Invitrogen)

Use a 1% DNA grade agarose (w/v) gel made with 0.5x TBE solution, and stained with 10% (v/v) SYBR safe.

6. 1x TE Buffer

| | Per 100 ml |
|-----------------------|------------|
| 1 M Tris-HCl (pH 8.0) | 1 ml |
| 0.5 M EDTA | 200 µl |

Adjust pH to 8.0± 0.2. Store at room temperature.

7. 6x loading dye

| | Per 100 ml |
|------------------|------------|
| 1 x TE | 10 ml |
| Glycerol | 50 ml |
| Bromophenol blue | 0.2% |

Store at room temperature.

Method

This method is to be repeated for each set of the three primer pairs listed in Table 1. Use one-step RT-PCR reagents as specified by the manufacturer. Some volumes outlined below may vary depending on the buffer and enzyme concentrations specified by the manufacturer. Ensure that the final volume of the RT-PCR reaction is 25 µl by altering the volume of nuclease-free water accordingly.

1. Label sterile PCR tubes
2. Prepare "Master Mix" on ice in a sterile microcentrifuge tube.

The "Master Mix" usually contains buffer, forward and reverse primers, RT/Taq and nuclease-free water.

Prepare the "Master Mix" according to the RT/Taq manufacturer's recommendations.

Ensure that the final volume for each reaction is 24 µl.

Add 24 µl of Master Mix to each PCR tube.

3. Add 1µl of template (total RNA extract) to the corresponding labelled PCR tube.
4. Cycle the tubes with the following PCR conditions: 1 cycle [48°C for 45 min], 1 cycle [94°C for 2 min], 35 cycles [92°C for 30 secs, annealing temperature as in Table 1 for 30 secs, 72°C for 60 sec], 1 cycle [72°C for 5 mins] and 15°C hold.

Please note that the PCR conditions have been adapted from Mayo et al. (1996) (25 cycles, annealing temperature of 50°C), Sokmen et al. (1998) (annealing temperature of 50°C), and Nassuth et al. (2000) (annealing temperature of 54°C for 45 seconds).

Cycling conditions for the nested-PCR round 1 are as follows: 1 cycle [48°C for 45 min], 1 cycle [94°C for 2 min], 40 cycles [92°C for 30 secs, annealing temperature as in 2 for 30 secs, 72°C for 60 sec], 1 cycle [72°C for 10 mins] and 15°C hold. Following completion of 1st round, follow manufacturer's instructions for DNA polymerase 2nd round PCR.

5. Cycling conditions for the nested-PCR round 2 are as follows: 1 cycle [94°C for 2 min], 40 cycles [92°C for 30 secs, annealing temperature as in 2 for 30 secs, 72°C for 60 sec], 1 cycle [72°C for 10 mins] and 15°C hold (Fig 19). At the completion of the RT-PCR, mix 10 µl each PCR sample with 2 µl of 6x gel loading dye, and load samples onto a 1% agarose gel with SYBR™ DNA Gel Stain (Invitrogen).

6. Electrophorese in 0.5X TBE at 100V for 45 minutes or until the bromophenol blue front has migrated half way down the length of the gel.
7. Visualise and photograph gel on UV transilluminator.
8. Criteria for determination of a valid assay: 196 bp product from samples containing plant extracts, no product in no template and negative plant extract controls, product of the expected size in positive control.

Criterion for positive sample result: Product of the expected size in test sample, obtained from a valid assay.

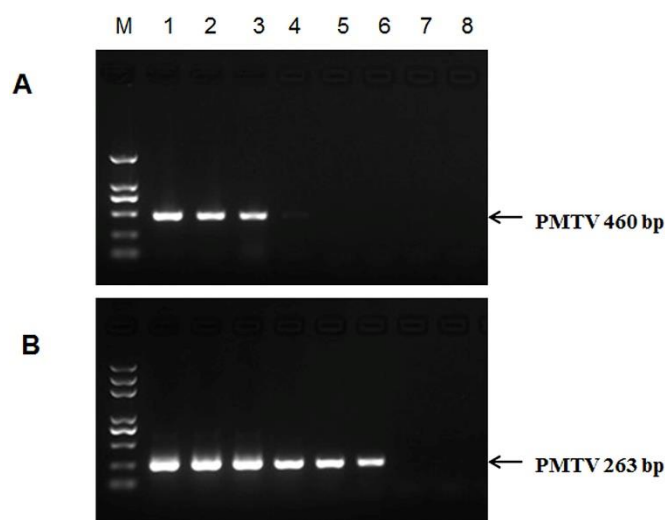


Fig. 19. The results of conventional RT-PCR and nested RT-PCR with serial RNA dilutions. Lane M, DNA marker; lanes 1–8 correspond to RNA dilutions of 100, 10⁻¹, 10⁻², 10⁻³, 10⁻⁴, 10⁻⁵, 10⁻⁶ and 10⁻⁷, respectively. A. Conventional RT-PCR. B. Nested RT-PCR using 0.5 μ L of the external PCR products as the template [Source: Zhou *et al.*, 2019].

4.1.3 DNA Sequencing of PCR Products

Sequencing is outsourced and the technique not described here. The following is the method used to prepare the sample for DNA sequencing.

Equipment required

1. 0-2 μ L, 2-20 μ L, 20-200 μ L, and 200-1000 μ L micropipettes and tips
2. 0.2 or 0.5 ml PCR tubes
3. 1.5 or 2 ml centrifuge tubes to store reagents
4. Bulb spinner or centrifuge
5. Freezer
6. Crushed ice
7. Latex gloves
8. PC with Internet access
9. Thermocycler
10. UV illuminator

Reagents

1. QIAquick PCR Purification Kit, or gel purification kit if multiple bands are present. Available from Qiagen, Catalogue Number 28104 (Alternatively any validated kit may be used). Forward and Reverse primer (As per section 0).
2. Sterile dH₂O.

Method

PCR products of correct size are purified using the QIAquick PCR Purification Kit (Qiagen) as per manufacturer's instructions. Sequencing can be out-sourced to a commercial provider (e.g. Micromon, Monash University; AGRF etc) and the DNA sequences are compared against sequences on the GenBank database (<http://www.ncbi.nlm.nih.gov>) using the program BlastN (Altschul *et al.*, 1997), to confirm if the positive PCR product sequence is from PMTV (PMTV RNA2 Genbank accession No. D30753; PMTV RNA 3 accession No. D16193).

4.2 Serological Tests – ELISA

The ELISA protocol for PMTV (Bioreba is the preferred manufacturer) detects all known strains of PMTV and does not cross react with other viruses.

For PMTV detection using double-antibody sandwich (DAS)-ELISA using the BIOREBA (Reinach, Switzerland) including interpretation of results, follow the manufacturer's instructions. For Further information: <http://www.bioreba.ch/saas/web/bioreba/web.aspx?PageID=58&search=PMTV>

The buffer recipes are include in Appendix 8.1.

4.2.1 Equipment

- 20, 100 and 1000ul pipettes and tips
- Microcentrifuge and microcentrifuge tubes (2.0mL)
- Container with ice
- Mortar and pestle (or similar macerating device)
- ELISA plates
- ELISA plate reader.

4.2.2 Reagents

- Coating Antibody (provided by supplier)
- Detecting Conjugate (provided by supplier)
- Controls: Positive and negative (provided by supplier – Permit required).
- 5X PBS Buffer
- PBS-Tween-Buffer
- Coating Buffer
- ELISA extraction Buffer
- Conjugation Buffer
- Substrate Buffer

5 CONTACTS FOR FURTHER INFORMATION

| AUSTRALIA | UNITED STATES OF AMERICA |
|--|--|
| <p>David Lovelock Department of Jobs, Precincts and Regions Agribio 5 Ring Road, La Trobe University Bundoora, Victoria 3083 Phone: 03 9032 7170 david.lovelock@agriculture.vic.gov.au</p> | <p>Steven B. Johnson Extension crops specialist Extension Professor University of Maine Cooperative Extension Aroostook City Ces Aroostook County Phone: 764-3361 stevenj@maine.edu</p> |
| CANADA | SCOTLAND |
| <p>Will Proctor Agriculture Information Officer 440 University Avenue Charlottetown Prince Edward Island C1A 4N6 Phone: (902) 620-3285 waproctor@gov.pe.ca</p> | <p>Lesley Torrance Scottish Crop Research Institute Invergowrie Dundee, DD2 5DA Scotland, UK Tel +44(0)1382 5627331 l.torrance@scri.sari.ac.uk</p> |

.....

6 ACKNOWLEDGEMENTS

This protocol was written by Jo Eyre, Bonny Rowles-van Rijswijk and Dr Brendan Rodoni from the Department of Primary Industries, Knoxfield, Private Mail Bag 15, Ferntree Gully Delivery Centre, Victoria 3156, Australia.

Dr John Thomas, AgriScience Queensland, peer reviewed and verified the protocol. It was reviewed and updated by Dr. David Lovelock at AgriBio.

7 REFERENCES

- Altschul SF, Madden TL, Sc  ffer AA, Zhang J, Zhang Z, Miller W and Lipman DJ (1997) Gapped BLAST and PSI-BLAST: a new generation of protein database search programs. *Nucleic Acids Research* **25**, 3389-3402.
- Andersen, BAB; Nicolaisen, M and Nielsen, SL (2002). Alternative hosts for *Potato mop-top virus*, genus *Pomovirus*, and its vector *Spongospora subterranea*. *Potato res* **45**, 37-43
- B  chen-Osmond C (2001 onwards) 27.0.1.0.004 *Potato mop-top virus*, In: ICTVdB, version 3, based on the 7th ICTV Report and subsequent up-dates.
<http://www.ictvdb.rothamsted.ac.uk/ICTVdB/27010004.htm>
- Calvert EL (1968) The reaction of potato varieties to *Potato mop-top virus*. *Res. exp. Rec. Minist. Agric. Nth. Ire.* **17**, 31-40.
- Cerovska N, Moravec T, Rosecka P, Dedic P, Filigarova M (2003) Production of polyclonal antibodies to a recombinant coat protein of *Potato mop-top virus*. *Journal of Phytopathology* **151**, 195-200.
- Cowan GH, Torrance L, Reavy B (1997) Detection of *Potato mop-top virus* capsid readthrough protein in virus particles. *Journal of General Virology* **78**, 1779-1783.
- Germundsson A, Sandgren M, Barker H, Savenkov EI, Valkonen JPT (2002) Initial infection of roots and leaves reveals different resistance phenotypes associated with coat protein gene-mediated resistance to *Potato mop-top virus*. *Journal of General Virology* **83**, 1201-1209.
- Gil, J.F., Adams, I., Boonham, N., Nielsen, S.L. and Nicolaisen, M. (2016), Molecular and biological characterization of Potato mop-top virus (PMTV, Pomovirus) isolates from the potato-growing regions of Colombia. *Plant Pathol*, 65: 1210-1220. doi:10.1111/ppa.12491.
- Harrison BD and Jones RAC (1970) Host range and some properties of *Potato mop-top virus*. *Annals of Applied Biology* **65**, 395-402.
- Helias V, Jacquot E, Guillet M, le Hingrat Y and Giblot-Ducray D (2003) Production of recombinant *Potato mop-top virus* coat protein in *Escherichia coli* and generation of antisera recognising native virus protein. *Journal of Virological Methods* **110**, 91-97.
- Johnson (2004) Potato Facts *Potato mop-top virus* (PMTV). *University of Maine Cooperative Extension Bulletin #2437* <http://www.umext.maine.edu/onlinepubs/htmlpubs/2437.htm>
- Jones RAC (1988) *Epidemiology and control of potato mop-top virus*. Dev. Appl. Biol. Vol. 2:255-270. **IN** Cooper JI and Asher MJC (editors) *Viruses with fungal vectors*. Proceedings of a conference at the University of St. Andrews, Aug. 25-27, 1987. Wellesbourne, Warwick, Great Britain.
- Jones RAC and Harrison BD (1969) The behaviour of *Potato mop-top virus* in soil and evidence for its transmission by *Spongospora subterranea* (Wallr.) Lagerh. *Annals of Applied Biology* **63**, 1-17.
- Jones RAC and Harrison BD (1972) Ecological studies on *Potato mop-top virus* in Scotland. *Annals of Applied Biology* **71**, 47-57.

- Kalyandurg, P., Gil, J.F., Lukhovitskaya, N.I., Flores, B., Müller, G., Chuquillanqui, C., Palomino, L., Monjane, A., Barker, I., Kreuze, J. and Savenkov, E.I. (2017), Molecular and pathobiological characterization of 61 Potato mop-top virus full-length cDNAs reveals great variability of the virus in the centre of potato domestication, novel genotypes and evidence for recombination. *Molecular Plant Pathology*, 18: 864-877. doi:10.1111/mpp.12552
- Kurppa AHJ (1989) Reaction of potato cultivars to primary and secondary infection by *Potato mop-top furovirus* and strategies for virus detection. *EPPO Bulletin* **19**, 593-598.
- MacKenzie DJ, McLean MA, Mukerji S, Green M (1997) Improved RNA extraction from woody plants for the detection of viral pathogens by reverse transcription - polymerase chain reaction. *Plant Disease* **81**, 222-226.
- Mayo MA, Torrance L, Cowan G, Jolly CA, Macintosh SM, Orrega R, Barrera C, Salazar LF (1996) Conservation of coat protein sequence among isolates of *Potato mop-top virus* from Scotland and Peru. *Archives of Virology* **141**, 1115-1121.
- Mumford RA, Walsh K, Barker I, Boonham N (2000) *Detection of Potato mop top virus and Tobacco rattle virus* using a multiplex real-time fluorescent reverse-transcription polymerase chain reaction assay. *Phytopathology* **90** (5), 448-453.
- Nassuth A, Pollari E, Helmezczy K, Stewart S and KoFalvi SA (2000) Improved RNA extraction and one-tube RT-PCR assay for simultaneous detection of control plant RNA plus several viruses in plant extracts. *Journal of Virological Methods* **90**:37-49.
- Nielsen SL, and Nicolaisen M. (2003) Identification of two nucleotide sequence sub-groups within *Potato mop-top virus*. *Archives of Virology* **148**, 381-388.
- Plant Health Risk Assessment Unit (2002) *Potato mop-top virus* (PMTV) <http://www.inspection.gc.ca/english/sci/surv/data/pmtvle.shtml>
- Santala, J., Samuilova, O., Hannukkala, A., Latvala, S., Kortemaa, H., Beuch, U., Kvarnheden, A., Persson, P., Topp, K., Ørstad, K., Spetz, C., Nielsen, S., Kirk, H., Budziszewska, M., Wieczorek, P., Obrępalska-Stęplowska, A., Pospieszny, H., Kryszczuk, A., Sztangret-Wiśniewska, J., Yin, Z., Chrzanowska, M., Zimnoch-Guzowska, E., Jackeviciene, E., Taluntytė, L., Pūpola, N., Mihailova, J., Lielmane, I., Järvekūlg, L., Kotkas, K., Rogozina, E., Sozonov, A., Tikhonovich, I., Horn, P., Broer, I., Kuusiene, S., Staniulis, J., Uth, J., Adam, G. and Valkonen, J. (2010), Detection, distribution and control of *Potato mop-top virus*, a soil-borne virus, in northern Europe. *Annals of Applied Biology*, 157: 163-178. doi:10.1111/j.1744-7348.2010.00423.x
- Scott KP, Kashiwazaki S, Reavy B, Harrison BD (1994) The nucleotide sequence of potato mop-top RNA 2: a novel type of genome organisation for a furovirus. *Journal of General Virology* **75**, 3561-3568.
- Sokmen MA, Barker H, and Torrance L (1998) Factors affecting the detection of *Potato mop-top virus* in potato tubers and improvement of test procedures for more reliable assays. *Annals of Applied Biology* **133**, 55-63.
- Tidona CA and Darai G (eds) (2002) *The Springer Index of Viruses*. Springer – Verlag Berlin, Germany, pg 1297-1301.

- Thompson JR, Wetzel S, Klerks MM et al., 2003. Multiplex RTPCR detection of four aphid-borne strawberry viruses in *Fragaria* spp. in combination with a plant mRNA specific internal control. *Journal of Virological Methods* 111, 85–93.
- Tsrer (Lahkim), L., Shapira, R., Erlich, O., Hazanovsky, M. and Lebiush, S. (2020), Characterization of weeds and rotational crops as alternative hosts of *Spongospora subterranea*, the causal agent of powdery scab in Israel. *Plant Pathol*, 69: 294-301. doi:10.1111/ppa.13117.
- Wicks T and de Boer D (2001) *Diseases – Issues and Control Practices*. South Australian Research and Development Institute, State of South Australia.
http://www.sardi.sa.gov.au/pages/horticulture/pathology/hort_pn_diseases.htm:sectID=334&tempID=111.
- Yellareddygari, S. K. R., Whitworth. L. J. & Gudmestad, N. C. (2018) Assessing Potato Cultivar Sensitivity to Tuber Necrosis Caused by Potato mop-top virus. *Plant Disease* **102**: 1148-1153.
- Zhou, H., Lei, Y., Wang, P., Liu, M. and Hu, X (2019) Development of SYBR Green real-time PCR and nested RT-PCR for the detection of *Potato Mop-top Virus* (PMTV) and viral surveys in Progeny tubers derived from PMTV infected Potato tubers. *Molecular and Cellular Probes* **47**, 101438.

7.1 Related Articles

- Arif M, Torrance L and Reavy B (1999) Acquisition and transmission of *Potato mop-top furovirus* by a culture of *Spongospora subterranea* f. sp. *subterranea* derived from a single cystosorus. *Annals of Applied Biology* **126**, 493-503.
- Calvert EL and Harrison BD (1966) Potato mop-top, a soil-borne virus. *Plant Pathology* **15**,134-139.
- Cooper JI and Harrison BD (1973) Distribution of *Potato mop-top virus* in Scotland in relation to soil and climate. *Plant Pathology* **22**(2), 73-78.
- Harrison BD and Jones RAC (1971) Factors affecting the development of spraing in potato tubers infected with *Potato mop-top virus*. *Annals of Applied Biology* **68**, 281-289.
- Hu X, Dickinson V, Changzheng, YL, Singh M, Yang Y, Xiong X and Nie X (2016) Molecular characterization of Potato mop-top virus isolates from China and Canada and development of RT-PCR differentiation of two sequence variant groups. *Canadian Journal of Plant Pathology* **38**, 231-242.
- Jones RAC and Harrison BD (1969) The behaviour of potato mop-top virus in soil and evidence for its transmission by *Spongospora subterranea* (Wallr.) Lagerh. *Annals of Applied Biology* **63**, 1-17.
- Kalischuk M, Lynn J and Kawchuk (2016) First report of Potato mop-top virus Infecting Potatoes in Alberta. *Plant Disease* 100:2544-2544. Doi:10.1094/PIDS-06-16-0902-PDN
- le Romancer M, Kerlan C and Nedellec M (1994) Biological characterisation of various geographical isolates of *Potato virus Y* inducing superficial necrosis on potato tubers. *Plant Pathology* **43**, 138-144.

- McGeachy KA, and Barker H (2000) *Potato mop-top virus* RNA can move long distance in the absence of coat protein: evidence from resistant, transgenic plants. *Molecular Plant-Microbe Interactions* **13** (1), 125-128.
- Reavy B, Sandgren M, Barker H, Heino P, Oxelfelt P (1997) A coat protein transgene from a Scottish isolate of *Potato mop-top virus* mediates strong resistance against Scandinavian isolates which have similar coat protein genes. *European Journal of Plant Pathology* **103** (9), 829-834.
- Sandgren M (1995) *Potato mop-top virus* (PMTV): distribution in Sweden, development of symptoms during storage and cultivar trials in field and glasshouse. *Potato Research* **38**, 379-389.
- Solomon-Blackburn RM and Barker H (2001) Breeding virus resistant potatoes (*Solanum tuberosum*): a review of traditional and molecular approaches. *Heredity* **86**, 17-35.
- Torrance L, Cowan GH, Arli Sokmen M and Reavy B (1999) A naturally occurring deleted form of RNA 2 of *Potato mop-top virus*. *Journal of General Virology* **80**, 2211-2215.

8 APPENDIX

8.1 ELISA Buffers

Coating buffer

Dissolve in distilled water to 1000 ml:

| | |
|------------------------------|--------|
| Sodium carbonate (anhydrous) | 1.59 g |
| Sodium hydrogen carbonate | 2.93 g |

Adjust pH to 9.6 and store at 4° C.

PBS-Tween Buffer (Wash Buffer)

Dissolve in distilled water to 1000 ml:

| | |
|--|--------|
| Sodium chloride | 8.0 g |
| Sodium phosphate, dibasic (anhydrous) | 1.15 g |
| Potassium phosphate, monobasic (anhydrous) | 0.2 g |
| Potassium chloride | 0.2 g |
| Tween-20 | 0.5 g |

Adjust pH to 7.4. Store at 4°C.

ECI Buffer (use as Sample Buffer and as Conjugate Buffer)

Add to 1000 ml 1X PBST:

| | |
|---|--------|
| Bovine serum albumin (BSA) | 2.0 g |
| Polyvinylpyrrolidone (PVP) MW 24-40,000 | 20.0 g |

Adjust pH to 7.4. Store at 4° C.

PNP Buffer

Dissolve in 800 ml distilled water:

| | |
|--------------------------------|---------|
| Magnesium chloride hexahydrate | 0.1 g |
| Sodium azide | 0.2 g |
| Diethanolamine | 97.0 ml |

Adjust pH to 9.8 with 0.1 N hydrochloric acid. Adjust final volume to 1000 ml with distilled water. Store at 4° C.

9 DIAGNOSTICS PROCEDURES TO SUPPORT SURVEILLANCE

9.1 Introduction

PMTV produces internal and external potato tuber symptoms, seriously affecting tuber quality and reducing yield (Germundsson *et al.*, 2002). Infection with PMTV causes some quantitative yield loss but the qualitative losses are more important and can lead to total crop rejection by supermarkets and processors (Jones and Harrison, 1972; Mumford *et al.*, 2000; Yellareddygaru *et al.*, 2018).

The current methods for detection include two RT-PCRs and ELISA and a newly added nested PCR. In situations where volumes of material/samples are very high (>300), it is recommended that the laboratory undertake ELISA initially, followed up with a more targeted approach using RT-PCR (Table 1).

Table 1: Diagnostic methods for the detection of PMTV.

| Method | Identification level | Identification Confidence | Deployment (Field/Lab) | Required Time | Throughput (No. of samples) |
|----------------|----------------------|---------------------------|------------------------|---------------|-----------------------------|
| RT-PCR (4.1.2) | To Species/Strain | High (99%+) | Laboratory | <1 w | High (100s) |
| ELISA (4.2) | To Species | Medium (90-99%) | Laboratory | <1 w | Very High (1000s) |

9.2 Sampling

Leaf and tuber samples should not have free water on their surfaces. Samples should not be collected during hot weather and all samples should be cooled and transported to a testing facility in a cool environment if possible. Further information can be found in the NDP sections 3.4.1 and 3.4.2.

9.3 In Field Tests

There are currently no known in field tests for PMTV.

9.4 Laboratory Tests

The laboratory tests including RT-PCR and ELISA can be found in the NDP for PMTV in sections 4.1.2 and 4.2 respectively.

9.5 Acknowledgements

This section was written and produced by David Lovelock, Agriculture Victoria and checked by the Diagnostic Protocol Working Group of SPHD.

9.6 References

- Germundsson A, Sandgren M, Barker H, Savenkov EI, Valkonen JPT (2002) Initial infection of roots and leaves reveals different resistance phenotypes associated with coat protein gene-mediated resistance to *Potato mop-top virus*. *Journal of General Virology* **83**, 1201-1209.
- Jones RAC and Harrison BD (1969) The behaviour of potato mop-top virus in soil and evidence for its transmission by *Spongospora subterranea* (Wallr.) Lagerh. *Annals of Applied Biology* **63**, 1-17.
- Mumford RA, Walsh K, Barker I, Boonham N (2000) *Detection of Potato mop top virus and Tobacco rattle virus* using a multiplex real-time fluorescent reverse-transcription polymerase chain reaction assay. *Phytopathology* **90** (5), 448-453.
- Yellareddygar, S. K. R., Whitworth. L. J. & Gudmestad, N. C. (2018) Assessing Potato Cultivar Sensitivity to Tuber Necrosis Caused by Potato mop-top virus. *Plant Disease* **102**: 1148-1153.